Safe
Heating and Mixing

Safety Guide Hei-PLATE

Index

German	3	
English	11	
French	18	
Español	26	
Italiano	34	
Português	42	
Română	50	
Magyar	58	
Hrvatski	65	
Български	72	
Ελληνική	80	
Slovenčina	88	
Slovenščina	95	
Polski	102	
Čeština	110	
Lietuvių kalba	117	
Eesti	124	
Latviešu	131	
Dansk	138	
Suomi	145	
Nederlands	152	
Svenska	160	

Zu diesem Dokument

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About this document

This document contains a summary of all safety instructions for Heidolph Scientific Products GmbH' magnetic stirrers and lab jacks.

This document in no way replaces the operating instructions supplied with the product!

Typographic conventions

Standardized symbols, signal words and highlighting elements are used in this document to warn of hazards and to identify important information or rather special text contents.

Symbol	Signal word / explanation
	Warning symbols in combination with a signal word indicate dangers:
	DANGER Indicates an immediate dangerous situation. Failure to respect the indications will result in death or serious injury.
	WARNING Indicates a potential danger. Failure to respect the indications will result in serious injuries.
	CAUTION Indicates a potential hazard which, If not avoided, damage to property and minor to moderate injuries can occur.
	Mandatory signs are used to indicate important and useful information on handling a product. This information is used to ensure operational safety and to maintain the value of the product.

Copyright protection

This document is protected by copyright and is intended for use by the purchaser of the product only.

No transfer to third parties, reproduction in any form, including excerpts, and by any means, as well as utilization and/or disclosure of the contents is permitted without the prior written consent of Heidolph Scientific Products GmbH. Any violation is subject to compensation for damage.

Intended use

Magnetic stirrers of the company Heidolph Scientific Products GmbH have been developed and designed for the following laboratory activities:

- Heating
- Mixing
- Stirring
- Titrating

Lab jacks of the company Heidolph Scientific Products GmbH have been exclusively developed and designed for the operation with magnetic stirrers, type Hei-PLATE Mix'n'Heat Ultimate (lifting and lowering of the magnetic stirrer placed on the lab jack).

Detailed information can be seen in the associated operating instructions. The areas of application of Heidolph Scientific Products GmbH' devices include especially chemical, biological and environmental-analytical laboratory and research applications, applications of basic research and comparable facilities.

Due to their design, the devices in their delivery conditions may only be used in analytical processes or in laboratory-like conditions in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals.

Any other use is deemed to be not as intended!

Compliant use

The user is generally responsible for evaluating the conformity of his application and, if necessary, for taking additional measures.

Reasonably foreseeable misuse

Additional measures may be necessary for use under conditions or for purposes deviating from the intended use, and/or specific guidelines and safety regulations must be observed. Corresponding requirements must be evaluated and implemented by the operator in each individual case.

Compliance with and implementation of all relevant directives and safety measures for the respective field of application is generally the responsibility of the operator.

All risks resulting from improper use are borne generally by the operator.

The device may generally be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the device is handled responsibly are the general responsibility of the operator!

Transportation

During transportation, avoid severe shocks and mechanical stresses that can cause damage to the device. Keep the original packaging in a dry and protected place for later use.

Storage

Always store the device in its original packaging. To protect against damage and excessive material aging, store the device in an environment that is as dry, temperature-stable and dust-free as possible.

Acclimatization

After each transportation and after storage under critical climatic conditions (e.g. high temperature difference between inside and outside), allow the device to acclimatize at room temperature for at least two hours to prevent possible damage due to condensation before commissioning it in the place of use. If necessary, extend the acclimatization phase if the temperature differences are very high.

Make all supply connections (power supply, tubing) only after the device has been acclimatized!

Permissible ambient conditions

The device is designed for indoor use only. The device is not suitable for outdoor use! The device is not suitable for use in potentially explosive areas!

When used in corrosive atmospheres, the service life of the device may be reduced depending on the concentration, duration and frequency of exposure.

General safety instructions

Before commissioning and using the device, familiarize yourself with all the safety regulations and occupational safety guidelines applicable at the place of use and observe them at all times.

Only operate the device if it is in faultless technical condition. In particular, ensure that there is no visible damage on the device itself and, where applicable, on connected devices or the supply connections.

If there is missing or misleading information on the device or regarding occupational safety, contact the responsible safety specialist or our technical service.

Only use the device in accordance with the regulations for intended use.

Electrical safety

Before connecting the device to the power supply, ensure that the voltage indicated on the rating plate matches the specifications of the local power utility company.

Ensure that the power supply circuit provided is protected by means of a residual-current device (RCD).

Always use the power supply cord supplied with the device. All installation work for the power supply of the device may only be carried out by an authorized electrician or by the technical service of Heidolph Scientific Products GmbH!

Prior to use, check that the device and the power supply cord are free of visible damage.

Always have repairs and/or maintenance work on the device carried out by an authorized electrician or by the technical service department of Heidolph Scientific Products GmbH.

Always switch off and disconnect the device from the power supply, preventing reconnection, before carrying out maintenance work, cleaning, or repairs.

Interfaces

Safely isolate extra low voltage inputs and outputs via 25 V AC or 60 V DC according to EN 61140, or by double or reinforced insulation according to EN 60730-1 or DIN 60950-1.

Use only shielded connection cables. Connect the shield to the connector housing.

Data security

The user is responsible for ensuring data security when transferring data between the described device and other devices.

Always use secure networks for the data transfer and avoid use of critical infrastructure.

Always use high-quality shielded data cables for the data transfer.

For data transfer via a USB B connection, an industrial standard USB hub should be preferably used to ensure the most stable connection possible.

Operational safety

Operate the device under a closed ventilated fume hood when working with potentially hazardous substances (see EN 14175 and DIN 12924).

Do not make any unauthorized changes or modifications to the device!

Always use genuine spare parts and accessories, or those expressly approved by the manufacturer!

Rectify malfunctions or faults on the device immediately.

Switch off and disconnect the device from the power supply, preventing reconnection, if it is not possible to eliminate the malfunction or rectify the fault immediately.

Observe all relevant general and safety instructions for the connected peripheral devices (observe the supplied documentation!).

Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.

Occupational safety

Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.

Persons with pacemakers or other susceptible implants must maintain a minimum distance of ten cm from the device to prevent possible interference.

Do not operate any other devices in the immediate vicinity of the device

- which generate emission or radiation sources in the frequency range 3×10^{11} Hz to 3×10^{15} Hz (in the optical spectral range wavelengths from 1,000 μm to 0,1 μm),
- which generate ultrasonic or ionizing waves.

Do not process any substances that could release energy in an uncontrolled manner (e.g. self-ignition).

Do not process substances in which the energy input through mixing poses a danger.

Do not work with an open flame in the immediate vicinity of the device (risk of explosion or fire when processing explosive/flammable media).

Wipe off any liquid that may have spilled on the device immediately.

Always switch off the device or disconnect the device from the power supply when it is not in use.

Personal protective equipment (PPE)

The operating company must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The corresponding instruction of the personnel is generally within the operating company's responsibility.

Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid hazards to the environment.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the general responsibility of the operator!

Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking the device with a biohazard warning symbol.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the general responsibility of the operator!

Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Noncompliance will invalidate any warranty claims against Heidolph Scientific Products GmbH.

The operator is generally liable for all damage resulting from unauthorized changes or modifications to the device, from the use of unapproved or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

Device-specific safety instructions (magnetic stirrer)

Mounting

CAUTION

To use the device, place it on a clean, stable, level, horizontal, and heat-resistant surface. Make sure that the power plug can be reached directly at any time, in order to disconnect the device from the power supply at any time without delay.

Before switching on, make sure that the device is sufficiently stable.

Before switching on, always make sure that all setups are centered as far as possible and are fixed on the base unit with sufficient stability.

Keep all support and contact surfaces clean and dry.

During operation, make sure that the required minimum distance of 15 cm between the magnetic stirrer and other devices and structures is maintained.

Power supply

ANGER

The device may only be supplied with the required mains voltage via a grounded mains socket-outlet (voltage see rating plate).

Only use the supplied power supply cord.

Operation

CAUTION

Whenever possible, use closed vessels for processing corrosive, toxic or biohazardous substances and seal them safely.

Increase the speed - especially at high speeds in combination with a large load and/or a tall setup on the heating plate - only gradually to the required level and pay attention to the stability of the setup.

Lower the speed or reduce the total load on the heating plate if the device starts to move independently during operation.

Always pay attention to even distribution of the containers and in particular to uniform distribution of the weight if vessels of different sizes and/or differently filled vessels are placed on the heating plate at the same time.

High pressures can result when processing samples in closed containers. Always use adequately dimensioned, pressure-resistant vessels: Note the expansion volume of the sample!

Always use the necessary and appropriate personal protective equipment!

Secure the device in remote mode with a clearly visible warning sign and, if necessary, take further ambient-specific protective measures that protect against damage to property and injuries in the event of unexpected/unmonitored starting of the device.

Heating mode

WARNING

Contact with surfaces above 50 °C can lead to serious injury.

Always pay attention to the heating temperature indicator of the device.

Avoid direct skin contact with the heating plate during operation and also observe the residual heat indicator after switching off!

Do not place heat-sensitive objects on the heating plate.

Always use the necessary personal protective equipment (heat-resistant gloves, eye protection, safety clothing) to process samples from a temperature of 50 °C.

As long as the device is not disconnected from the power supply, the display shows the actual temperature of the heating plate or temperature measured at the external temperature sensor after switching off the device (residual heat indicator).

Note that the temperature at the external sensor can differ substantially from the surface temperature of the heating plate.

Avoid direct skin contact with the heating plate!

Do not place heat-sensitive objects on the heating plate.

Rotational mode

WARNING

When using open vessels, there is a risk that fluid will spray out.

Whenever possible, use closed vessels for processing corrosive, toxic or biohazardous substances and seal them safely.

Especially using open vessels, increase the rotation speed only gradually and observe the fluid movements.

Always use the necessary personal protective equipment (heat-resistant gloves, eye protection, safety clothing) to process samples from a temperature of 50 °C.

Automatic restart

WARNING

When the [automatic restart] function is activated, the heating plate automatically heats up to the last set operating temperature when the power is restored and the rotation speed will be augmented to the last set value.

Ensure that all vessels are properly positioned on the heating plate before the power is restored. In the event of a power failure, do not place any objects on the heating plate.

Magnetic coupling interrupted detection function

CAUTION

In the event of a magnetic coupling interruption, there is a risk of the magnetic stirring bar thrashing about in the container in an uncontrolled way and damaging the glass.

Only deactivate the magnetic coupling interrupted detection function if this is useful for the process or is required.

Operation with external temperature sensor/temperature control unit

WARNING

Note that if a temperature sensor is connected, the temperature is adjusted to it automatically. As long as the temperature sensor is not immersed in the sample, the reference temperature of the ambient air is displayed. The heating plate can therefore heat up to 350 °C unnoticed!

Always immerse the temperature sensor in the sample before switching on the device or before calibrating!

Do not place heat-sensitive objects on the heating plate.

Always use the necessary personal protective equipment (heat-resistant gloves, eye protection, safety clothing) to process samples from a temperature of 50 °C.

Remote mode

WARNING

Secure the device in remote mode with a clearly visible warning sign and, if necessary, take further ambient-specific protective measures that protect against damage to property and injuries in the event of unexpected/unmonitored starting of the device.

Device service

DANGER

Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

When cleaning, avoid the penetration of liquids.

Before replacing the fuses, switch off the device and disconnect the power supply cord.

Always replace the two device fuses in pairs with original manufacturer fuses. Further information on available accessories can be found on our website at www.heidolph.com!

After fuse replacement, check the device for a safe condition according to IEC 61010-1.

CAUTION

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

Device-specific safety instructions (lab jack)

CAUTION

The USB interface type C of the magnetic stirrer may only be used for operation with a Heidolph Scientific Products GmbH' lab jack. The general USB-C standard is not supported. Direct or indirect damage caused by connecting devices not authorized by the manufacturer is the responsibility of the user.

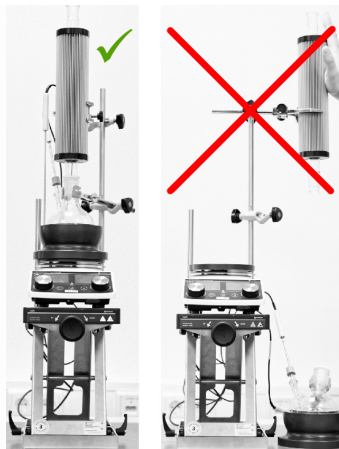
WARNING

The lab jack is designed for a load of maximum 10 kg. Direct or indirect damages caused by an overload of the lab jack, are generally the sole responsibility of the user.

Always be sure to position all setups and accessory parts centrally above the magnetic stirrer on the lab jack.

Move the lab jack downward and/or pull the setups centrally upward as much as possible to remove glass flasks or attachments from the magnetic stirrer!

Never pivot setups to the outside (see example: setup with magnetic stirrer, heating attachment and Findenser®).

Mounting

CAUTION

Observe the instructions for the correct positioning of the device.

Observe the generally valid rules for the safe installation of work equipment.

To use the lab jack, place it on a clean, dry, stable, even and horizontal surface.

Always make sure that all setup components and accessory parts of the magnetic stirrer are securely fixed (see also the corresponding delivery documentation or operating and/or mounting instructions):

Fix the setups preferably on a wall grid (if available). When mounted on a wall grid, ensure that the wall grid and all mounting elements have sufficient self-protection and load capacity!

Use all recommended retaining and fixing elements for setup components and accessory parts.

The lab jack, the magnetic stirrer and any other setups do not form a physical unit and must always be transported separately.

Before each transport, disconnect the cable connection between the lab jack and the magnetic stirrer.

The professional mounting and positioning of the device including the supplied accessories is within the sole responsibility of the operator!

Heidolph Scientific Products GmbH accepts no responsibility for direct and/or indirect personal injury or damage to property resulting from failure to observe the instructions for the correct mounting and positioning of the device.

Operation

CAUTION

Never reach into the lab jack mechanism after it is connected to the magnetic stirrer.

Do not place any objects in the lab jack mechanism.

Always make sure that there is sufficient distance between the lab jack and other devices or structures: the required minimum distance is 15 cm on all sides.

Make sure that no incoming or outgoing lines (power connection cable, data cable, hose connections) protrude into the lab jack's mechanism.

Clean the lab jack's mechanism at regular intervals or immediately if it is highly soiled.

Always make sure that the travel path of the lab jack is not blocked.

In the event of high setups, pay particular attention to moving the center of gravity of the whole setup to the middle of the lab jack table as far as possible to ensure adequate stability.

Attach a suitable warning of unexpected traversing movements of the lab jack to the overall setup in remote and in unsupervised operation.

Device service

CAUTION

In the event of contact with aggressive media, make sure that all surfaces are cleaned immediately with suitable aids to prevent damage to the device surface.

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

Repairs

Repairs to the device may only be carried out by authorized experts!

Unauthorized repairs during the warranty period will result in the loss of the warranty claim.

The owner is solely liable for damage caused by unauthorized repairs.

If repairs are necessary, contact an authorized dealer or our technical service.

Maintenance

The device contains no user-serviceable components. If necessary (in the event of abnormal operating behavior such as excessive noise or heat generation, for example) contact our technical service.

Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Contact details

Heidolph Scientific Products GmbH

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© Heidolph Scientific Products GmbH