Safety Guide

Hei-VAP

Index

Deutsch	2	
English	11	
Български	20	
Čeština	30	
Tysk	39	
ελληνικά	47	
Español	58	
Eesti	67	
Suomi	76	
Français	85	
Hrvatski	94	
Magyar	103	
Italiano	112	
Lietuvių k.....	121	
Latviešu	130	
Nederlands	139	
Polski	148	
Português	157	
Română	166	
Slovenčina	175	
slovensko	184	
Svenska	194	

Zu diesem Dokument

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www.heidolph.com

About this document

This document contains a summary of all safety instructions for the following devices:

- Hei-VOLUME Distimatic Pro
- Hei-VAP Expert/Ultimate
- Hei-VAP Industrial/HBX

This document in no way replaces the operating instructions supplied with the product! For more detailed information, please refer to the original instructions (in German) or the translation of the original instructions supplied with the product.

Typographic conventions

The following symbols, signal words and highlighting elements are used in this document:

Symbol	Signal word / explanation
	Warning symbols in combination with a signal word indicate dangers: DANGER Indicates an immediate dangerous situation. Failure to respect the indications will result in death or serious injury. WARNING Indicates a potential danger. Failure to respect the indications will result in serious injuries. CAUTION Indicates a potential hazard which, if not avoided, damage to property and minor to moderate injuries can occur.
	Mandatory signs are used to indicate important and useful information on handling a product. This information is used to ensure operational safety and to maintain the value of the product.

Copyright protection

This document is protected by copyright and is generally intended for use by the purchaser of the product.

No transfer to third parties, reproduction in any form, including excerpts, and by any means, as well as utilization and/or disclosure of the contents is permitted without the prior written consent of Heidolph Scientific Products GmbH. Any violation is subject to compensation for damage.

Customs declaration

Rotary evaporators, as combustion and cleaning equipment, may be subject to notification to a competent customs authority in the country of destination.

The evaluation of a customer obligation to notify and, if applicable, the notification to a competent customs authority in the country of destination is generally the responsibility of the user!

Intended use

The described device was designed by the manufacturer for the separation of substances, powder drying, concentration, crystallization of substances, and recycling of solvents under vacuum.

Due to its design, the device in its delivery condition may generally be used in analytical processes or in laboratory-like conditions in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals.

Any other use of this device is not considered as intended!

Compliant use

The user is generally responsible for evaluating the conformity of his application and, if necessary, for taking additional measures.

Reasonably foreseeable misuse

Additional measures may be necessary for use under conditions or for purposes deviating from the intended use, and/or specific guidelines and safety regulations must be observed. Corresponding requirements must be evaluated and implemented by the operator in each individual case.

Compliance with and implementation of all relevant directives and safety measures for the respective field of application is generally the responsibility of the operator.

All risks resulting from improper use are borne generally by the operator.

The device may generally be operated by authorized and instructed personnel. Training and qualification of the operating personnel as well as ensuring that the device is handled responsibly are the general responsibility of the operator!

Transportation

During transportation, avoid severe shocks and mechanical stresses that can cause damage to the device. Keep the original packaging in a dry and protected place for later use.

Storage

Always store the device in its original packaging. To protect against damage and excessive material aging, store the device in an environment that is as dry, temperature-stable and dust-free as possible.

Acclimatization

After each transportation and after storage under critical climatic conditions (e.g. high temperature difference between inside and outside), allow the device to acclimatize at room temperature for at least two hours to prevent possible damage due to condensation before commissioning it in the place of use. If necessary, extend the acclimatization phase if the temperature differences are very high.

Make all supply connections (power supply, tubing) only after the device has been acclimatized!

Permissible ambient conditions

The device is designed for indoor use only. The device is not suitable for outdoor use! The device is not suitable for use in potentially explosive areas!

When used in corrosive atmospheres, the service life of the device may be reduced depending on the concentration, duration and frequency of exposure.

General safety instructions

Before commissioning and using the device, familiarize yourself with all the safety regulations and occupational safety guidelines applicable at the place of use and observe them at all times.

Only operate the device if it is in faultless technical condition. In particular, ensure that there is no visible damage on the device itself and, where applicable, on connected devices or the supply connections.

If there is missing or misleading information on the device or regarding occupational safety, contact the responsible safety specialist or our technical service.

Only use the device in accordance with the regulations for intended use.

Electrical safety

Before connecting the device to the power supply, ensure that the voltage indicated on the rating plate matches the specifications of the local power utility company.

Ensure that the power supply circuit provided is protected by means of a residual-current device (RCD).

Always use the power supply cord supplied with the device. All installation work for the power supply of the device may only be carried out by an authorized electrician or by the technical service of Heidolph Scientific Products GmbH!

Prior to use, check that the device and the power supply cord are free of visible damage.

Always have repairs and/or maintenance work on the device carried out by an authorized electrician or by the technical service department of Heidolph Scientific Products GmbH.

Always switch off and disconnect the device from the power supply, preventing reconnection, before carrying out maintenance work, cleaning, or repairs.

Data security

The user is responsible for ensuring data security when transferring data between the described device and other devices.

Always use secure networks for the data transfer and avoid use of critical infrastructure.

Always use high-quality shielded data cables for the data transfer.

For data transfer via a USB B connection, an industrial standard USB hub should be preferably used to ensure the most stable connection possible.

Operational safety

Operate the device under a closed ventilated fume hood when working with potentially hazardous substances (see EN 14175 and DIN 12924).

Do not make any unauthorized changes or modifications to the device!

Always use genuine spare parts and accessories, or those expressly approved by the manufacturer!

Rectify malfunctions or faults on the device immediately.

Switch off and disconnect the device from the power supply, preventing reconnection, if it is not possible to eliminate the malfunction or rectify the fault immediately.

Observe all relevant general and safety instructions for the connected peripheral devices (observe the supplied documentation!).

Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.

Occupational safety

Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.

Do not operate any other devices in the immediate vicinity of the device ...

- which can generate electromagnetic fields in the frequency range between 9×10^3 Hz to 3×10^{11} Hz,
- which generate emission or radiation sources in the frequency range 3×10^{11} Hz to 3×10^{15} Hz (in the optical spectral range wavelengths from 1,000 μm to 0,1 μm),
- which generate ultrasonic or ionizing waves.

Do not operate the unit when adiabatic compression or shock waves may occur (pressure wave ignition).

Do not use substances that could release energy in an uncontrolled way and cause a pressure increase (exothermic reaction, spontaneous ignition of dusts).

Do not process hard, brittle materials such as stones, soil samples, etc., that could destroy the evaporation flask.

Only use heating bath media that guarantee sufficient heat transfer.

Do not operate the device with overpressure.

Do not expose the glass components to a pressure difference of more than 2 bar.

Ensure that the coolant overpressure does not exceed a level of 2 bar.

Ensure that the flow velocity does not exceed 1 m/s when aspirating liquids with flammable components (electrostatic charge, danger of ignition!).

Avoid the formation of explosion group IIC gases and potentially explosive distillation residues.

Personal protective equipment (PPE)

The operating company must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The corresponding instruction of the personnel is generally within the operating company's responsibility.

Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid hazards to the environment.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the general responsibility of the operator!

Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking the device with a biohazard warning symbol.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the relevant personnel is the general responsibility of the operator!

Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

Noncompliance will invalidate any warranty claims against Heidolph Scientific Products GmbH.

The operator is generally liable for all damage resulting from unauthorized changes or modifications to the device, from the use of unapproved or non-genuine spare parts and accessories, or from disregarding the safety instructions and hazard warnings or the manufacturer's instructions!

Specific safety instructions Hei-VOLUME Distimatic Pro

Assembly

CAUTION: Material damage, loss of production

Due to improper system tubing and wiring as well as unauthorized changes to the system, there is a risk of direct and indirect damage to property!

For system tubing and wiring, follow all notes and instructions provided in the assembly instructions provided (part number 01-001-009-12-0).

Do not make any unauthorized changes to the device, system, tubing, nor the existing cable connections.

If necessary, contact our technical service (contact details see operating instructions).

Control panel

CAUTION: Damage to the control panel

Tip or sharp objects can damage the surface of the display on the control panel!

To operate the device, touch the display only with your fingertips or a special stylus with a soft-touch tip.

Switch the device on/off

CAUTION: Property damage, loss of production

If the control box is switched on and off incorrectly, there is a risk that peripheral devices will not be detected correctly or will switch in an uncontrolled manner!

Observe the order given when switching the control box and the connected peripheral devices on and off.

In particular, before turning off the control box, make sure that all processes have been completed properly.

Observe the following sequence when turning the automatic module and attached peripherals on and off:

TURNING ON THE SYSTEM: First, turn on all peripheral devices so that they are recognized when initializing the automatic module software.

TURNING OFF THE SYSTEM: First, turn off all peripherals to put the entire system in a safe operating condition.

Calibrate the flask level sensor

CAUTION: Loss of production

The function of the level sensor in the evaporation flask (S4) may be affected by electromagnetic fields. This can lead to a loss of the operational quality of the entire system.

During calibration, commissioning and operation, ensure that there are no sources of interference in the immediate vicinity of the evaporation flask to ensure proper operation of the level sensor at all times.

System check

CAUTION: Property damage, loss of production

Faulty connections or signaling can lead to malfunctions or uncontrolled switching operations!

Leaks reduce the performance of the entire system.

Correct any indicated errors and/or detected malfunctions immediately.

In the event of a continued fault indication/malfunction, contact your local sales representative or our technical service department (contact details see operating instructions).

Perform overpressure test

WARNING: Risk Of Injury, Damage To Property

Due to faulty components, faulty assembly of components or faulty connections, there is a risk that pressure-applied components will suddenly loosen or rupture during an overpressure test!

Before performing an overpressure test, ensure that all components are properly installed.

In particular, make sure that all screw connections are properly tightened.

Before starting the overpressure test, close all safety covers, hoods and doors of the entire system.

Ensure that third parties are clear of the danger zone.

Use the required personal protective equipment (eye protection, hand protection, etc.).

Specific safety instructions Hei-VAP Industrial

Control panel

CAUTION

Pointed or sharp objects can damage the surface of the display on the control panel!

To operate the device functions, always touch the display with your fingertips or a special stylus with a soft-touch tip.

Interfaces

WARNING

If the interfaces are assigned too high voltages and are not sufficiently insulated, metallic parts such as the housing can be live in the event of a fault.

Safely separate low-voltage inputs and outputs via 25 V AC or 60 V DC according to DIN EN 61140, or by double or reinforced insulation according to DIN EN 60730-1 or DIN 60950-1.

Always use shielded connection cables. Connect the shield to the connector housing.

CAUTION

Make sure that the device interfaces on the bottom of the display are always properly covered!

Penetrating liquid can cause interface malfunctions and short-circuiting.

Cooling water supply

CAUTION

Keep the [Supply] and [Return condenser] connections closed if no external condenser is connected!

In order to not exceed the maximum permissible pressure in the glass condenser (2 bar) during operation, open the [Supply] and [Return condenser] connections when an external condenser is connected!

EASY LOCK flask clamping system

CAUTION

Incorrect or careless handling can damage the flask clamping system and/or the flask inserted!

Before inserting a flask for the first time, familiarize yourself with the handling of the EASY LOCK flask clamping system.

Before inserting a flask for the first time, familiarize yourself with the handling of the flask support.

Unsuitable flasks can damage the flask clamping system and/or the flask inserted!

The EASY LOCK flask clamping system is only suitable for flasks with an opening diameter of 150 mm. The use of flasks with a smaller opening diameter (e.g. LR 20) is not permitted, as these cannot be fixed in a positive manner due to the geometric properties.

Excessive contact/closing pressure can damage the EASY LOCK and/or the flask inserted!

Always follow the instructions for the correct adjustment of the contact pressure.

Heating bath

DANGER

When using oils with a flash point < 285 °C as a heating bath liquid, uncontrolled thermal reactions can occur at high heating bath temperatures and an explosive atmosphere can arise: Always use oils with a flash point > 285 °C as heating bath liquid.

From a heating bath temperature of approx. 75 °C – 80 °C, the use of polyethylene glycol or oil is recommended as heating bath liquid.

CAUTION

When the evaporation flask is immersed, the already filled heating bath liquid is displaced and the liquid level in the heating bath rises!

Especially when using oil, the device and the surrounding floor surface can be heavily contaminated by overflowing heating bath liquid.

When filling the heating bath, make sure that the maximum filling quantity is not exceeded.

The monitoring of the refill duration during the automatic water bath refill should only be deactivated in exceptional cases. (only Hei-VAP Industrial)

Consider the mark for maximum filling level in the heating bath or the information for maximum filling volume when choosing the manual refill!

CAUTION

The heating bath is designed for a maximum operating temperature of 210°C (using silicone oil). From a liquid temperature of 50 °C, there is a risk of injury in the event of contact.

Always grip and hold the heating bath by the ergonomic safety handles on the side! (only Hei-VAP Expert/Ultimate)

Before emptying the heating bath, allow heated liquid to cool down to below 50 °C.

Use suitable personal protective equipment (heat-resistant gloves, eye protection, safety clothing) to empty the heating bath.

Guard hood (optional)

DANGER

Even when the guard hood is closed, there is a risk that the heating bath liquid will spill out of the heating bath at high rotational speeds. There is a risk of scalding and/or contamination.

CAUTION

Always use the handle provided to open and close the guard hood.

When closing the guard hood, keep sufficient distance to the closing edges with your free hand.

Set up the device

CAUTION

Due to improper mounting and/or installation of the system/device, as well as unauthorized changes to the mounting, there is a risk of direct and indirect damage to property!

If necessary (system mounting and installation carried out by the operator or third parties, necessary changes to the existing mounting), contact our technical service.

To use the device, place it on a clean, stable, level and horizontal surface.

Before switching on the device, make sure that it is sufficiently stable and use all available locking means on the base.

Specific safety instructions Hei-VAP Expert/Ultimate

Control panel

CAUTION

Pointed or sharp objects can damage the surface of the display on the control panel!

To operate the device functions, always touch the display with your fingertips or a special stylus with a soft-touch tip.

Interfaces

CAUTION

Penetrating liquid can cause malfunctions and short-circuiting.
Open the cap for the lateral interfaces only if needed and make sure that it is always connected in normal operation.

Set up the device

CAUTION

Due to improper mounting and/or installation of the system/device, as well as unauthorized changes to the mounting, there is a risk of direct and indirect damage to property! If necessary (system mounting and installation carried out by the operator or third parties, necessary changes to the existing mounting), contact our technical service.
To use the device, place it on a clean, stable, level and horizontal surface.

PTFE seal

CAUTION

Improper cleaning can damage the surfaces of the PTFE seal.
Clean the seal surfaces with a soft, lint-free and only slightly moistened cloth.
Never use any aggressive or abrasive cleaning agents or aids.

Safety instructions for rotary evaporators in general

Switching the device on/off

CAUTION

Before each operation of the device, check the caps and connecting elements of all glass components for correct and tight fit!
At loose/leaky connections, ambient air is sucked in during the vacuum build-up and the required vacuum pressure cannot be reached in a stable manner!
In the event of excess pressure, there is a risk that liquid will escape from the system! The device and surrounding areas can be severely contaminated by leaking media!
Tighten loose caps/fasteners by hand or correct the seating of the retaining clips.

Mounting/removing evaporation flasks

WARNING

In the event of improper handling, the evaporation flask that is to be used may fall to the ground.

The smallest damage to the evaporation flask can lead to glass breakage when pressurized.

- Glass breakage!
- Risk of injury/contamination by leaking medium!

Use only the evaporation flasks approved by the manufacturer.

When inserting and removing an evaporation flask, observe all instructions for the correct handling of the respective support system.

Before each operation of the device, check the evaporation flask for visible damage.

Operating modes

WARNING

Due to incorrect/insufficient process settings, there is a risk that impermissible operating values are reached and, as a result, device components will be damaged and heating bath liquid and/or distillation material will escape.

In general, qualified personnel is allowed to adjust the process settings of the device.

When determining the pressure values, observe the chemical-physical properties of the respective material to be distilled.

Pay particular attention to the specific safety instructions for the processing of toxic substances!

When operating the device, there is a risk of contact with rotating parts. Loose clothing, jewelry and open hair can be pulled in!

If possible, operate the device with a guard hood (accessory) or under a closed laboratory fume hood.

Never touch rotating parts with your hands. Do not bend over rotating parts.

Wear tight-fitting work clothes, remove jewelry before starting work, tie long hair under a hair net or under a suitable head covering.

Use appropriate personal protective equipment (PPE).

At high rotation speeds, the heating bath liquid can spray away from the evaporation flask by forming film and/or slop out of the heating bath. There is a risk of scalding and/or contamination.

If possible, operate the device with the guard hood closed.

Adjust the rotation speed of the evaporation flask according to the recommendations in this section.

Use appropriate personal protective equipment (PPE).

Clean dirty surfaces immediately!

Device parameters

CAUTION

Due to incorrect/insufficient safety settings there is a risk that processes will not be shut down or will be shut down with a time delay when an unacceptable pressure level is reached.

Only qualified personnel may adjust the safety settings of the device.

If necessary, contact our technical service.

Distillation under atmospheric pressure

CAUTION

During distillation under atmospheric pressure, excess pressure may occur in the system. There is a risk that, in a high system pressure situation, gas or liquid will escape into the ambient air in the area of the seals and joints. Risk of personal injuries/damage to property

There is a risk of bursting under extreme pressure conditions!

Carefully compare the process parameters with the prevailing permissible ambient conditions and the chemical-physical properties of the respective material to be distilled (refer to data sheets)!

Pay particular attention to the specific safety instructions for the processing of toxic substances!

Only operate the device with the guard hood closed or under a closed laboratory fume hood.

Use appropriate personal protective equipment (PPE).

Remote mode

WARNING

Secure the device in remote mode with a clearly visible warning sign and, if necessary, take further ambient-specific protective measures that protect against damage to property and injuries in the event of unexpected/unmonitored starting of the device.

Set rotation speed

WARNING

When operating the device without a guard hood, there is a risk of contact with rotating parts. Loose clothing, jewelry and open hair can be pulled in!

If possible, operate the device with the guard hood closed or under a closed laboratory fume hood.

Never touch rotating parts with your hands. Do not bend over rotating parts.

Wear tight-fitting work clothes, remove jewelry before starting work, tie long hair under a hair net or under a suitable head covering.

At high rotation speeds, the heating bath liquid can spray away from the evaporation flask by forming film and/or slop out of the heating bath. There is a risk of scalding and/or contamination.

If possible, operate the device with the guard hood closed or under a closed laboratory fume hood.

Adjust the rotation speed of the evaporation flask according to the manufacturer's recommendations.

Use appropriate personal protective equipment (PPE).

Clean dirty surfaces immediately!

Even when the guard hood is closed, there is a risk that the heating bath liquid will spill out of the heating bath at high rotational speeds. There is a risk of scalding and/or contamination.

Adjust the rotation speed of the evaporation flask according to the manufacturer's recommendations.

Use appropriate personal protective equipment (PPE).

Clean dirty surfaces immediately!

Device service

DANGER

Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

When cleaning, avoid the penetration of liquids.

Before replacing the fuses, switch off the device and disconnect the power supply.

Always replace the two device fuses in pairs with original manufacturer fuses. Further information on available accessories can be found on our website at www.heidolph.com!

After fuse replacement, check the device for a safe condition according to IEC 61010-1.

CAUTION

Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

Never use any aggressive or abrasive cleaning agents or aids.

Repairs

Repairs to the device may always be carried out by authorized experts! Unauthorized repairs during the warranty period will result in the loss of the warranty claim. The owner is generally liable for damage caused by unauthorized repairs. If repairs are necessary, contact an authorized dealer or our technical service.

Maintenance

There are no user-serviceable components in the unit housing. If necessary (in the event of abnormal operating behavior such as excessive noise or heat generation, for example) contact our technical service.

Check the caps and fasteners of all glass components for correct and tight fit at regular intervals and especially before each start-up of the device.

Check the level sensor in the heating bath regularly for proper operation, contamination, and limescale build-up to prevent malfunctions.

CAUTION

At loose/leaky connections, ambient air is sucked in during the vacuum build-up and the required vacuum pressure cannot be reached in a stable manner!

In the event of excess pressure, there is a risk that liquid will escape from the system! The device and surrounding areas can be severely contaminated by leaking media!

Tighten loose caps/fasteners by hand or correct the seating of the retaining clips.

Dismantling

Observe the instructions given in the associated assembly instructions when dismantling the device.

Disposal

When disposing of the device, observe the provisions of the WEEE Directive 2012/19/EU and its transposition into national law in the country of use.

When disposing of portable batteries, observe the provisions of the European Battery Directive 2013/56/EU and their implementation in national law in the country of use.

Check the device and all components for residues of substances that are hazardous to health, the environment and biohazardous before disposal.

Properly remove and dispose of residues of substances that are hazardous to health, the environment and biohazardous!

Contact details

Heidolph Scientific Products GmbH

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© Heidolph Instruments GmbH & Co. KG